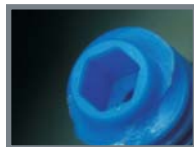
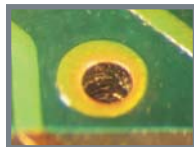
NEU ++ NEW ++ NEU
Mantis Elite-Cam
Imaging Software

Vision
ENGINEERING

Vifox Imaging

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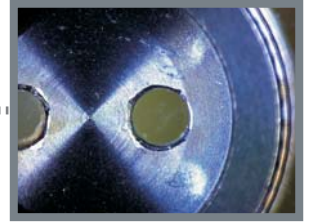
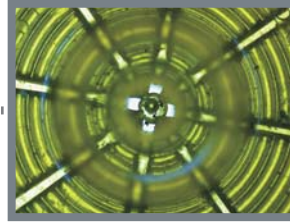
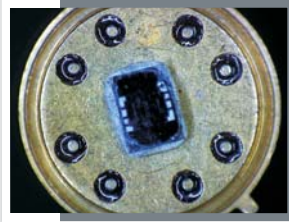
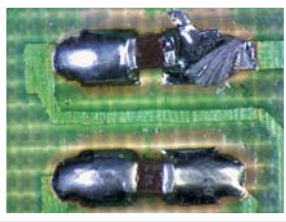
*Vifox is an
imaging software package for Mantis Elite,
allowing easy image acquisition,
measurement and storage of images.*

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Vifox is an image acquisition and measurement software package easily integrated with the Mantis Elite-Cam stereo viewer, providing precise imaging of your samples and components.

A single cable conveniently handles the data transfer - no separate mains adapter or additional wires!

The intuitive Vifox software allows you to easily set key parameters whilst viewing the live image. Then simply capture, document and archive your images.

The best pictures in the shortest time!

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Features:

- + On-screen settings for immediate results.
- + Images can be saved as *.jpg, *.bmp, *.png
- + Crosshair displayed as overlay
- + Software for measuring (2D) of lines, circles, areas, angles, drop perpendicular and more.
- + Scalebar in various displays, online value preview, digital zoom, show ruler, colour settings for lines and labelling, draw arrows and text, layer functionality
- + Stage micrometer for calibration purposes

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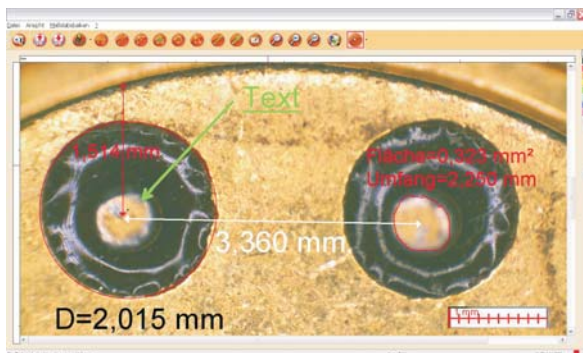
*You would like to measure image contents two-dimensional (2D)?
No problem with the measuring module of Vifox...*


1. Calibrate the system with the provided stage micrometer (< 1 minute).
2. Measuring unit and decimal places can be set during calibration.
3. Settings are stored on the PC for future use.
4. Acquire an image and assign the appropriate calibration.
- 5...The system is now ready to measure! (e.g. lines, circles, areas, angles, ...)

*Straightforward and easy to use, ideal for daily routine applications.
User-friendly features such as on-screen measuring value preview and digital zoom, allow precise setting of measurement points. Additional features include, scale bars, arrows and labelling, with the innovative layering function making further editing easy!*

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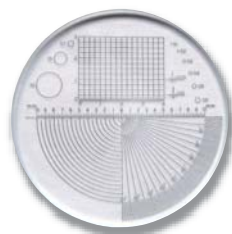*2D measurement function:
Comprehensive measuring options,
clear graphic presentation,
simple calibration, with
scale bars and labelling.*

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*Stage micrometer for calibration
purposes (included in delivery)*

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All function and operation steps in Vifox for measuring, camera settings and image acquisition are documented in a detailed HTML based online help section (F1 key). Comprehensive and illustrated installation instructions in PDF format are included on the product CD-ROM.

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Would you like to find out more about Vifox?

Visit our website at www.visioneng.com/vifox
...or ask your relevant sales consultant or distributor.

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Vifox wurde entwickelt von unserem Imaging-Partner
dhs Solution GmbH | www.dhssolution.com
Vifox was developed by our imaging partner
dhs Solution GmbH | www.dhssolution.com

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Vifox is available in English, German, French, Italian and Spanish languages and sold exclusively by Vision Engineering and authorised distributors.

Vision Engineering and authorised Distributors can provide technical support for Vifox software.

For Vifox system requirements, please visit www.visioneng.com/vifox

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von Vision Engineering Ltd.
Vision Engineering Ltd.